

The Picus Pipette Quick Guide

PICUS PIPETTE SETUP PROCEDURE

Ten Volumes Can Be Programed On The Sartorius Picus Pipette (10-300uL)

1. Use the **ON / OFF** Button to turn on the Pipette **(B)**
2. Press the **TIP EJECT** Button to pre-set to Ready **(D)**
3. Using the **SOFTKEYS**, press **MENU**, then **PIPETTE**, then **OK** **(F)**
4. Press **EDIT** using the **SOFTKEYS** (*Volume Will Be Highlighted*) **(F)**
5. Use the Black **ADJUSTMENT WHEEL** around Orange Operating Button to set new volume **(C)**
6. Press **NEXT** using the **MIDDLE SOFTKEY** to select speed (*Directions set speed at 5↑, 6↓*) **(F)**
7. Press **OK** once speed is selected using the **MIDDLE SOFTKEY** **(F)**
8. Press the **HOTKEY** Button above the Softkey Menu Buttons to select saved location **(E)**
9. **Scroll** using the black **ADJUSTMENT WHEEL** to desired location **(C)**
10. Using the **SOFTKEYS**, press **SAVE** and **EDIT** to name the location **(F)**
11. Use the black **ADJUSTMENT WHEEL** **(C)** to select chosen letter or press the **MIDDLE SOFTKEY** **(F)** to advance to Numbers / Clear
12. Press the Orange **OPERATING BUTTON** to advance to next programming position **(A)**
13. Press **OK** using the **MIDDLE SOFTKEY** when finished **(F)**



RECOMMENDED PROGRAMABLE VOLUMES

Program Positions	M1	M2	M3	M4	M5	M6
Volume uL	225	25	200	250	150	300