
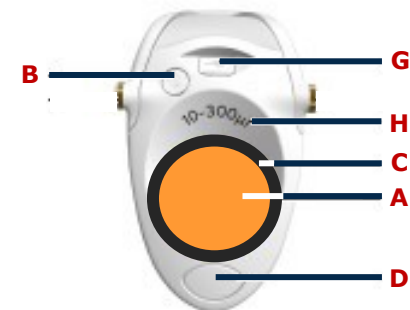


# The PICUS<sup>®</sup> 2 Pipette Quick Guide

## Programmable Volume and Memory Set Up Procedure

1. Use the **ON / OFF (B)** Button to Turn ON the Pipette
2. Press the **TIP EJECT (D)** Button
3. Press the **RIGHT SOFTKEY (F)** to Display Application Menu
4. PIPETTING Should be Highlighted, if Not Use the Black **SCROLLWHEEL (C)** to Select Pipetting
5. Press OK Using the **RIGHT SOFTKEY (F)**
6. Press EDIT Using the **MIDDLE SOFTKEY (F)** to Set Volume
7. Use the Black **SCROLLWHEEL (C)** to Select New Volume
8. Press NEXT Using the **MIDDLE SOFTKEY (F)** to Set Aspirate Speed
9. Use the Black **SCROLLWHEEL (C)** to Set Aspirate Speed to 5 ^
10. Press NEXT Using the **MIDDLE SOFTKEY (F)** to Set Dispense Speed
11. Use the Black **SCROLLWHEEL (C)** to Set Dispense Speed to 6 v
12. Press OK Once Speeds are Entered Using the **RIGHT SOFTKEY (F)**
13. Press the **HOTKEY (E)** to View Memory Locations
14. Use the Black **SCROLLWHEEL (C)** to Select the Desired Memory Location
15. Press the **MIDDLE SOFTKEY (F)** to SAVE
16. Using the **MIDDLE SOFTKEY (F)**, Press EDIT to Name the Memory Location
17. Use the Black **SCROLLWHEEL (C)** to Select Letters to Name Memory Location  
(Press the **MIDDLE SOFTKEY (F)** to Switch to Lowercase Letters, Numbers, Symbols or to Clear Entry)
18. To Save Each Letter or Number Selected, Press OK using the **RIGHT SOFTKEY (F)** to Move to the Next Space
19. After Entering the Name, Press OK Repeatedly Using the **RIGHT SOFTKEY (F)** Until the Curser Disappears
20. Press SAVE Using the **RIGHT SOFTKEY (F)**; the Entered Name is Now Stored in the Memory
21. Press the Orange **OPERATING BUTTON (A)** to Begin Pipetting Under this Memory Location
22. Repeat the Steps Above to Program Up to 20 Memory Locations

RECOMMENDED PROGRAMABLE VOLUMES & MEMORY LOCATIONS						
Memory Location	M1	M2	M3	M4	M5	M6
Volume (µL)	225	25	200	250	150	300

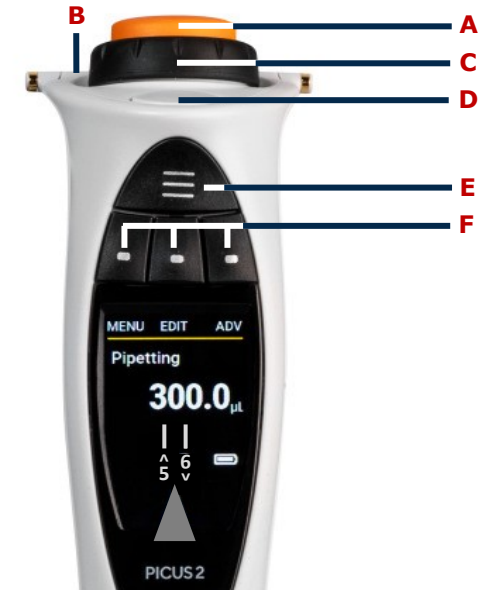

**NOTE:** The **CABLE CHARGING PORT (G)** and the **PIPETTE VOLUME RANGE (H)** locations are also detailed. PICUS<sup>®</sup> is a Registered Trademark of Sartorius Lab Instruments GmbH.

155 Gibraltar Road  
Horsham, PA 19044 USA

**Worldwide:** +1 215-441-4000  
**USA:** 1-800-257-3282  
**Fax Worldwide:** +1 215-443-8820  
**customer.service@biodatacorp.com**

**REF**  
**106075**  
**106077**  
**106232**

 **BIO/DATA**  
**CORPORATION**  
An ISO 13485 Registered Company  
[www.biodatacorp.com](http://www.biodatacorp.com)

Proudly Manufactured in the USA

**EC** **REP**

mdi Europa GmbH  
Langenhagener Str. 71  
D-30855 Langenhagen GERMANY

**UK** **REP**

Alpha Laboratories  
40 Parham Drive Eastleigh  
SO50 4NU Hampshire UNITED KINGDOM

